

Midazolam and 1-Hydroxymidazolam in Human Plasma : LDTD-MS/MS Analysis in 8 seconds

Pierre Picard and Patrice Tremblay

Keywords: High-throughput, LDTD, Tandem mass spectrometry, Midazolam, 1-Hydroxymidazolam, Human Plasma

Overview

- High-throughput determination of midazolam and 1-hydroxymidazolam in human plasma is performed by LDTD-MS/MS;
- Calibration range from 0.5 to 250 ng/mL with $r^2 > 0.997$;
- Sample-to-sample run time of 8 seconds;
- Accuracy within 94.9 and 112 %;
- Precision within 1.3 and 10.9 %.

Instrumentation

- Phytronix Technologies LDTD ion source (model T-960);
- Thermo Fisher Scientific TSQ[®] Quantum[™] Ultra AM mass spectrometer.

LDTD ionization process

The LDTD ion source uses an infrared laser diode to desorb sample that have been dried onto a well of a LazWell[™] (96-well plate). The desorbed gas phase molecules are carried into a corona discharge region to undergo APCI, then they are transferred directly into the mass spectrometer for detection.

Samples Preparation

Human plasma was spiked with midazolam and 1-hydroxymidazolam and with ISTD (deuterated midazolam and 1-hydroxymidazolam). The drugs were extracted with MTBE and hexane (3:1 v/v) and reconstitute into a water:acetonitrile solution (75:25 v/v). A volume of 2.0 μ L was manually transferred into a well of a LazWell[™] and was allowed to dry at room temperature.

Results and Discussion

Calibration Curves

Quantitative determination of midazolam and 1-hydroxymidazolam in human plasma can be achieved over a nominal concentration range of 0.5 to 250 ng/mL (**Figure 1 and 2**). An excellent linearity is obtained over the concentration range ($R^2 > 0.997$).

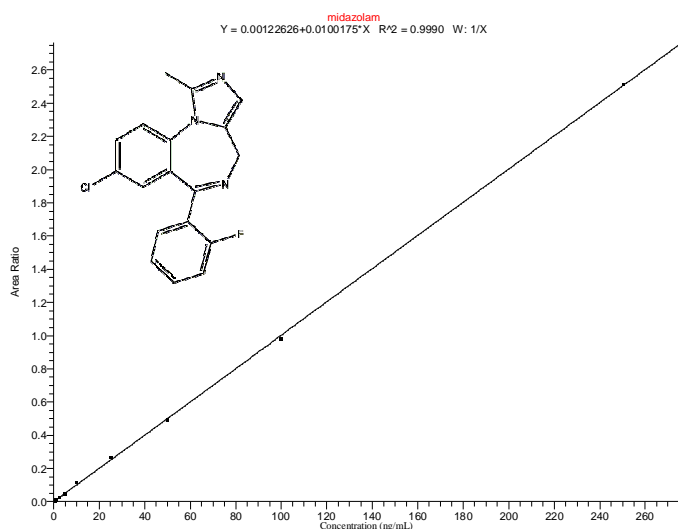


Figure 1 Calibration curve of midazolam in human plasma.

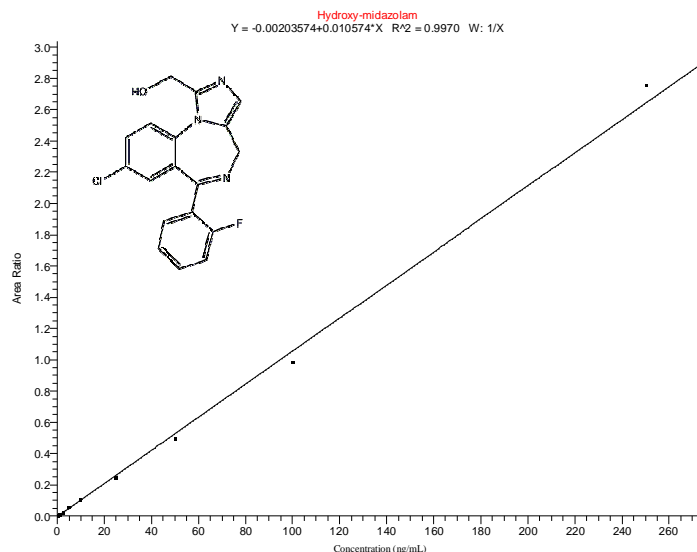


Figure 2 Calibration curve of 1-hydroxymidazolam in human plasma.

Accuracy and Precision

Quality control samples were analyzed in triplicates to evaluate the LDTD-MS/MS method accuracy and precision. The accuracy was evaluated to be within 94.9 and 112.2 % and the precision was within 1.3 and 10.9 % (**Table 1**)

Table 1 Within-run accuracy and precision for midazolam and 1-hydroxymidazolam.

	Midazolam		
	QC1	QC2	QC3
Nominal conc. (ng/mL)	1.5	15	200
N	3	3	3
Mean (ng/mL)	1.42	14.3	190.0
RSD (%)	10.9	6.0	1.3
% Nominal conc.	94.9	95.0	95.0
	1-Hydroxymidazolam		
	QC1	QC2	QC3
Nominal conc. (ng/mL)	1.5	15	200
N	3	3	3
Mean (ng/mL)	1.68	15.7	191.6
RSD (%)	3.8	4.3	1.6
% Nominal conc.	112.2	104.8	95.8

Midazolam and 1-Hydroxymidazolam LDTD Desorption Profile

The LDTD allows fast midazolam and 1-hydroxymidazolam thermal desorption (**Figure 3**). Blank sample signals allow a LOQ of 0.1 ng/mL for midazolam and of 0.8 ng/mL for 1-hydroxymidazolam.

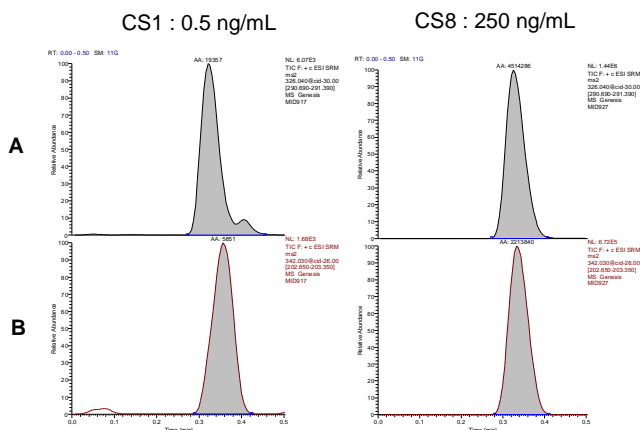


Figure 3 Desorption profile for 2 calibration standards for A) midazolam, and B) 1-hydroxymidazolam.

1-Hydroxymidazolam Back-conversion Evaluation

1-Hydroxymidazolam may back-convert into midazolam during the thermal desorption and/or the ionization process. To evaluate this back-conversion, 5 ng of 1-hydroxymidazolam was analyzed in LDTD-MS/MS and both midazolam and 1-hydroxymidazolam SRM transitions were monitored. The back-conversion of 1-hydroxymidazolam into midazolam was evaluated to be 0.18 % which does not affect midazolam quantification at plasma concentration.

MS Parameters

APCI (+)	
Collision gas pressure	1.5 mTorr (Argon)
Collision energy	27 V
Tube lens	77 V
Scan time	0.050 s
Needle voltage	5000 V
Q1 width	0.70 amu
Q3 width	0.70 amu
Midazolam SRM transition	326.04 → 291.04 amu
Midazolam-d4 SRM transition	330.00 → 295.00 amu
1-OH-midazolam SRM transition	342.03 → 203.00 amu
1-OH-midazolam-d4 SRM transition	346.00 → 203.00 amu

LDTD Parameters

Laser power pattern	0 to 40 % in 2.0 s Hold at 40 % for 2.0 s
Carrier gas flow	3.0 L/min (Air)

Conclusions

LDTD-MS/MS allows high-throughput quantification of midazolam and 1-hydroxymidazolam with a sample-to-sample run time of 8 seconds. Low back-conversion (0.8 %) was observed on 1-hydroxymidazolam.

High-throughput analysis with excellent linearity, accuracy and precision can be achieved using LDTD as ion source in mass spectrometry.

For more information about your specific application, visit www.phytronix.com

Phytronix
Technologies

337 rue Saint-Joseph E.
Québec, Qc Canada G1K 3B3
Phone : +1 (418) 692-1414
Fax : +1 (418) 692-4940

Copyright © 2008
All rights reserved